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Measure up

Thank you for purchasing a **separate ORP measurement electrode** or a **separate argentometry measurement electrode**.

For the best results with your electrode:

- **read** this user manual carefully,
- **comply** with the precautions for use.

The ORP electrode is used to measure a solution's oxidation reduction potential.
ORP = Oxidation Reduction Potential.

The separate argentometry electrode is used in association with a reference electrode to measure the concentration of silver ions in a solution.

The CE marking indicates compliance with the European directives, particularly LV and EMC.

The crossed-out dustbin means that, in the European Union, the product is the subject of selective disposal as per directive DEEE 2002/96/EC: this equipment must not be treated as household waste.

USE

- Remove the protective cap.
- Connect the reference electrode and measurement electrode to the instrument.
- Open the opening tab on the reference electrode.
- Calibration is not necessary. To test your electrode, use ORP test solutions.
- Immerse the tips of the electrodes in the solution to be measured, to the same height, while keeping them vertical and making sure that the membrane of the measurement electrode and the diaphragm of the reference electrode are immersed, but without exceeding the level of the internal electrolyte.
- Wait for the measurement to stabilize.
- Rinse the electrode with demineralized water after each measurement.
- Close the tab on the reference electrode again when you have finished measuring.

CLEANING

Metal electrodes may be soiled by absorption or oxidation. Clean them with demineralized water every time before and after use and then wipe them off with a clean, lint-free cloth.

If necessary, complete by washing with a dilute acid (such as acetic acid or HCl 0.01 mol/L) or a solvent (ethanol).

In some cases, it may be necessary to use a slightly abrasive cloth to remove persistent deposits. However, great care should be taken when using this method to avoid damaging the electrode.

Do not use aggressive products (hydrofluoric acid, acetone, etc.).

MAINTENANCE

Regularly check that your electrode remains clean and in good condition.

Keep your electrode in dry conditions with its protective cap.

Electrode reference	Filling electrolyte	Storage solution
BPT1, XPT1/XPT2	None	In dry conditions with protective cap
BAG1, XAG1		

STATE AT DELIVERY

The separate ORP measurement electrode or separate argentometry electrode is delivered in a cardboard box with:

- a shim,
- a protective cap,
- a bilingual user's manual,
- a multilingual accompanying datasheet.

For accessories and replacement parts, please visit our website:

www.chauvin-arnoux.com

LIFE SPAN AND WARRANTY

Separate ORP measurement electrodes and separate argentometry electrodes are wear parts. The life span of your electrode depends on the conditions in which it is used and the maintenance which you carry out on it.

The warranty does not apply if the glass is broken.

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